

General Tips and Tricks

Native MS

BEN JONES AND SOPHIE HARVEY
January 2022

Keep in mind the following tips and tricks from the Wysocki group when performing native mass spectrometry experiments.

#1: Tip Size Matters

- Different samples will require different kinds of tips for optimal nanospray.
- Larger tips which use higher spray voltage (1-1.4 kV) often work better for larger systems.
- Small (<150 kDa) soluble globular proteins tend to spray best from small tips at low voltage (0.5-0.9 kV).
- Tip size and spray voltage will drastically affect signal and apparent resolution or desolvation
- Don't be afraid to have multiple tip profiles.

#2: Use Fresh Solution of Cesium Iodide

- Cesium iodide for high mass calibration does not keep in solution. Prepare fresh for each calibration in 50/50 water/isopropanol at 2-8 mg/mL dependent on the instrument
- We recommend dissolving ~100 mg of CsI in a small amount of water (~1 mL), aliquoting (20 uL for 2 mg) into 1.5 - 2 mL tubes and then drying for storage. When needed you can resuspend an aliquot.

#3: Blunt Tipped Syringe for Filling Nanospray Capillaries

- If filling Nanospray capillaries using a syringe a blunt tipped syringe is recommended (Hamilton 80085 or similar).

#4: Consider Gel-Loading Tips

- Gel-loading tips (Genesse # 14-352, Eppendorf cat # 022351656) are useful for samples that may contaminate a syringe (ligands, detergent, salt, etc) and to eliminate the possibility of carryover.

continued...

#5: Recovering Limited Sample

- Sample can be recovered from a used tip, dispensed into a small tube, and reloaded into a new tip if sample is extremely limited.
- Samples that are sensitive to air or have been sprayed at high voltage (>2 kV) should not be transferred due to potential oxidation or other degradation.
- Attempting to directly transfer solution tip to tip often results in air bubbles that disrupt spray.

#6: Frequent Source Cleaning

- Simple source cleaning (not requiring venting) should be done daily for best results-particularly when additives such as magnesium acetate or detergents are being used. Frequent source cleaning can reduce internal contamination.
- Some samples are best run using a separate cone or ion transfer tube if available e.g. membrane proteins in detergent.
- Clean Pt wire often, or if contamination is suspected.

#7: Preferred Electrolyte Solutions

- Native mass spec electrolyte solutions are commonly ammonium acetate (Sigma 431311), ethylene diamine diacetate (EDDA) (Sigma 420352) [*Note: Not Ethylenediamine-N,N'-diacetic acid*], and triethyl ammonium acetate (TEAA) (Sigma 90358)
- Commonly used at 100 and 200 mM concentrations
- Complex formation may be affected by ionic strength and pH
 - Screening conditions may be required for unknown samples
- TEAA is added to 20% ionic strength in ammonium acetate ex 160 mM AmAc + 40 mM TEAA to achieve charge reduction
- Recommended to dilute a small amount of 1 M TEAA stock with water to typical ammonium acetate concentration for easy addition of 20% V/V to achieve charge reduction

#8: Using Ammonium acetate

- Ammonium acetate is hygroscopic and quickly absorbs water from the air making accurate ionic strength determination difficult. We recommend dissolving the entire container when received to 5 M in ultrapure water, storing at -20 C, and diluting immediately before use
- Ammonium acetate is volatile and concentration will change over time if left at room temperature or uncapped for extended periods of time
- It is recommended to dilute fresh ammonium acetate every 2 weeks for best results

#9: Dilute, Dilute, Dilute!

- Protein or protein complex concentrations should generally be 500 nM – 10 uM for analysis. Poor spray stability can be an indication of too high concentration, diluting 2-20x is recommended for testing. Excessive concentration can also result in the formation of nonspecific oligomers that are not present in solution